**INORA**  
Pharmaceutical Machinery

**INORA**  
Pharmaceutical Machinery

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# 01

Introduction

**INORA**  
Pharmaceutical Machinery Expert

Inora is a unique professional machine manufacturer in Taiwan. We engaged in researching, developing, designing, manufacturing, assembling, and consulting services for biotechnology and pharmaceutical industry. Our products and services are widely used in the pharmaceutical, food, biotechnology, chemical, and cosmetic fields which include mill granulator equipment, mixer, injection line, ointment line, extraction and concentration equipment. The precision and quality of its products have been recognized by all users.

Inora will keep servicing with the best management philosophy which suggests another landmark and the successful expansion to the global market of the company.

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## QUALITY

At Inora, quality inspection and assurance are an essential parts of the manufacturing process. We will take every possible measure to keep our product quality and customer satisfaction, so you don't have to worry about it, because we really understands what quality means to you.

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### THEORY +

Inside the cone-shaped vessel, the main impeller mix the powder into fluidized vortex and form lumps with the binder solution. then the high speed chopper cut the lumps into even granules by high speed action.

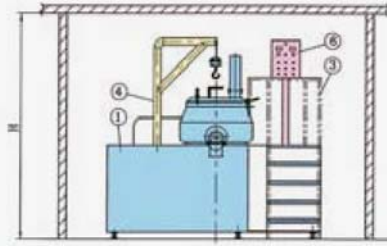
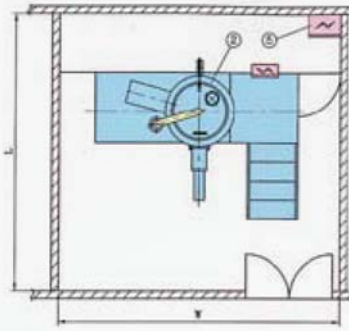
### APPLICATION +

- **Bio-tech / Pharmaceutical:** powder mixing and granulating before tableting.
- **Chinese medicine:** powder and soakage plaster mixing.
- **Food:** seasoning mixing and granulating.
- **Chemical:**
  - Chemical material mixing and granulating.
  - Plastic powder mixing and granulating.
  - Electronic metal powder mixing and granulating.

### FUNCTION +

- **Mixing:** powder and powder.
- **Granulation:** powder and binder solution.

## 流程圖 | FLOWCHART +

- 1. Frame
- 2. Mixer Tank
- 3. Ladder
- 4. Elevator
- 5. Control system
- 6. Operation Box

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## FEATURE +

- High uniformity, very different proportion also can uniformly mixing.
- Short mixing time, about 5 ~ 10 min/batch finished.
- Good tightness, use air bag to clamp the cover.
- Multi safety interlock device, high safety.
- Optional equipment for different product feature.
- Multi functions of mixing, granulating and drying can produce hormone or anti-cancer medicine.
- All the design and manufacture are qualified cGMP standard and FDA requirement.

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## THEORY +

Inside the cone-shaped vessel, the main impeller mix the powder into fluidized vortex and form lumps with the binder solution. then the high speed chopper cut the lumps into even granules by high speed action.

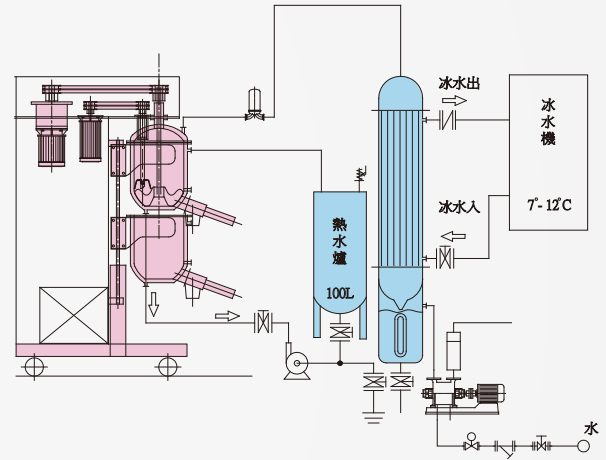
## APPLICATION +

- **Bio-tech / Pharmaceutical:**  
powder mixing and granulating before tableting.
- **Chinese medicine:**  
powder and soakage plaster mixing.
- **Food:**  
seasoning mixing and granulating.
- **Chemical:**
  - Chemical material mixing and granulating.
  - Electronic metal powder mixing and granulating.
  - Plastic powder mixing and granulating.

## FEATURES +

- Explosion Proof System is available: Suitable for organic solvent
- WIP System is available
- Multi safety interlock device, high safety.
- The clearance between impeller and bottom is very thin, therefore no dead angles and less remainder.
- Unique air seal system, a filtered compressed air is introduced into the air seal system. It prevents the powder fall into shaft and chopper, which may cause black spot and pollute the product.
- Good tightness, use air packing to clamp the cover..
- Multi functions of mixing, granulating and drying can produce hormone or anti-cancer medicine.
- All the design and manufacture are qualified cGMP , PIC/S standard and FDA requirement.

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## AUTOMATIC/CLOSED TYPE OPERATION +

- Powder Vacuum Loading is available
- Binder Spraying: Using the mist nozzle design
- Touch Screen Interface control system: all the parameters can be memorable, multilevel of pins can be set, connect to the diagrams printing

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Is used mainly in mixed of powder

### FEATURE +

- No screw in the mixer for greatly reducing pollution.
- Reversible running combined with double-layer adverse flow. Fast mixing. Convenient material infeed.

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### FUNCTION +

Granulate the wet powder into long granule.

### APPLICATION +

Pharmaceutical, food, chemical.

### FEATURE +

- Produced in a large number, fine granule, the size of granule uniformly.
- Screen diameter can be changed to adjust granule size.
- Making granule continuously.
- Convenient and precise for cleaning.
- All the design and manufacture are qualified cGMP standard and FDA requirement.

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## THEORY +

Two categories of granulation methods : **WET & DRY**

### WET TYPE:

granulation requires water or solvent to be added during the granulation process; consequently a drying process is essential after granulation stage.

### DRY TYPE:

granulation eliminates the water, or solvent adding process, and as a result the drying process is not required. However binder or lubricant can still be added during the drying granulation process.

Some raw material have the characteristic of low density (low specific gravity) and bad flow ability. By using the Roller Compactor-dry granulation process, it will improve the above characteristics and result in a bigger bulk density, higher specific gravity and good flow ability of granules.

## FEATURE +

- Three functions in one unit compression, milling and granulating.
- No water or solvent is added, granules are highly stable after granulation.
- Without water adding in the process, drying process is not required. Therefore, granulation process is easier.
- The good flow ability of outcome granules can be easily filled into capsules.
- Option to equip a vibration sifter, it filtrates small or undersize granules for re-granulation purpose.
- An automatic suction conveyer for charging powder or granule is available as an option for saving labor cost.

## APPLICATION +

- **Pharmaceutical:** Chinese Medicine 、 Western Drug 、 Antibiotic Chemical 、 Agriculture Product 、 Food.
- Producing granules of higher specific gravity and better flow ability.

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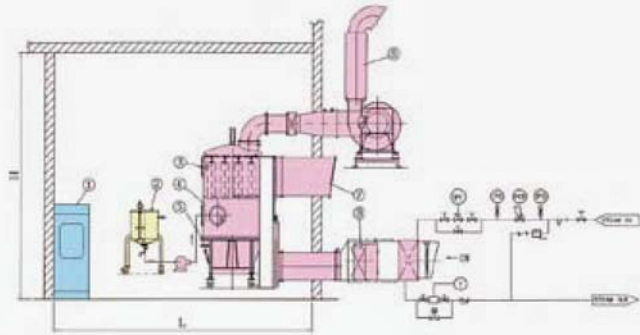
### THEORY +

Heated and purified air flows into the column by suction fan, the pellet or powder will be in uniform fluid condition inside the column, use one set of filter bags to prevent leak of powder, discharge the evaporated water to get draying purpose. If add one more spray gun can spray the binder or active ingredient downward into the uniform fluid powder, whenever the powder combined in granule will be dried to get the required moisture content.

### APPLICATION +

- **Bio-tech / Pharmaceutical:**
  - A. Western medicine: granulating and drying before tableting.
  - B. Chinese medicine: granulating and drying of concentrated granule.
- **Food:** Granulating and drying of instant food, instant health food, coffee, cocoa, milk powder, fruit flavor powder, seasoning.
- **Chemical:** Granulating and drying of pigment, pigment, dye, synthesis raw material.

- 1. Control Panel
- 2. Solution Tank
- 3. Filter
- 4. Drying Body
- 5. Trolley
- 6. Centrifugal Fan
- 7. Exhaust pipe
- 8. Heater

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## FEATURE +

- Mixing, granulating, drying are in one machine.
- Drying time of wet granule is 10 ~ 20 times faster than conventional oven.
- No dead angel produce because of even fluid drying, no color changing of difference because of the same drying speed in any position.

## FUNCTION +

- **Drying**: drying for powder or granule which contains moisture.
- **Granulating / Drying**: spray binder or active ingredients on tiny powder to combine in granules, drying until required moisture content.

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型號 Model	台車全容積 (L)	實際操作量 (KG) 依假比重:0.3計算	外型尺寸 Size (MM) (W/L/H)
YS-FDG-2	7	0.5-02	2500X2500X2100
YS-FDG-5	26	02-05	2500X3000X2300
YS-FDG-60	180	20-60	4000X4000X2800
YS-FDG-120	400	50-120	4000X4000X3200
YS-FDG-200	650	60-200	4500X5000X4500
YS-FDG-400	1300	100-400	5000X5000X5000
YS-FDG-600	2000	200-600	6000X6000X6000

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
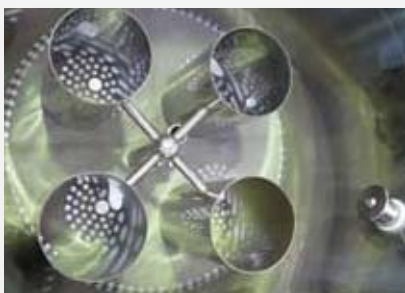
### FUNCTION +

- Top spray type:  
Drying wet granule or powder; granulation and drying of spraying binder on tiny powder.
- Bottom spray type :  
Spray suspension on pellet or film / enteric coating.
- Side spray type:  
Pellet powder coating; suspension coating or film / enteric coating.

### APPLICATION +

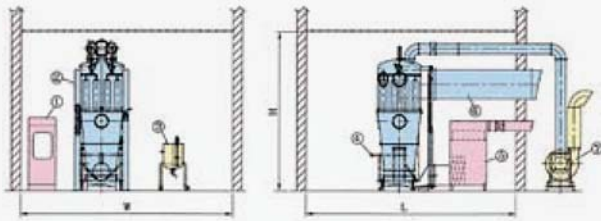- Top spray type:
  - Pharmaceutical: granulating and drying of Chinese or Western medicine, Bio-tech products.
  - Instant food: granulating and drying for instant powder of health food; juice extract powder; meat extract powder; seasoning; lactic acid bacteria.
  - Chemical: drying of dye and synthesis raw material.
- Bottom & Side spray type:  
Produce pellet of medicine and health food.

- 1. Control Panel
- 2. Drying Body
- 3. Solution Tank
- 4. Trolley
- 5. Heater
- 6. Exhaust pipe
- 7. Centrifugal Fan

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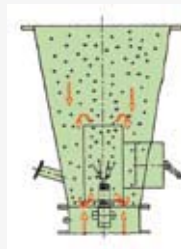
頂噴式  
Top spray type


discharge the vaporated water to get drying purpose. If add one more spray gun can spray the binder or active ingredient downward into the uniform fluid powder, whenever the powder combined in granule, will be dried to get the required moisture content.

- **Bottom spray type:**  
When the hot air flows through the bottom screen of coating column, it will generate siphonal principle. Convection is created through the strong force the granule will be sucked into the coating column and floating from bottom toward top, then fall down again, the bottom spray gun will spray upward into the uniform fluid granule to get coating and drying purpose.

底噴式  
Bottom spray type

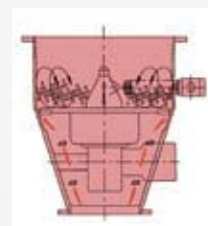- **Side spray type:**  
Place the sugar cores on the turntable. Hot air is blown upward between turntable and granulation area. The air causes the cores to roll, the binder solution is sprayed on the rolling cores through the pump and spray gun, coating and drying simultaneously. If powder coating is required, the fixed volume powder will be sprayed at the same time let solution coating the powder on the cores to produce pellet and get coating and drying purpose. cores to roll, the binder solution is sprayed on the rolling cores through the pump and spray gun, coating and drying simultaneously.

側噴式  
Side spray type

If powder coating is required, the fixed volume powder will be sprayed at the same time let solution coating the powder on the cores to produce pellet and get coating and drying purpose.

## THEORY +

- **Top spray type:**  
Heated and purified air flows into the column by suction fan, the pellet or powder will be in uniform fluid condition inside the column, use one set of filter bag to prevent leak of powder,

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型號 Model	台車全容積 Full Trolley Volume (L)	實際操作量 (KG) 依假比重:0.3計算	外型尺寸 Size (MM) (W/L/H)
YS-FD-2BG	7	0.5-2.0	2500X2500X2600
YS-FD-5BG	26	02-05	2500X3000X2600
YS-FD-60BG	180	20-60	4000X4000X3500
YS-FD-120BG	400	50-120	4000X4000X4000
YS-FD-200BG	650	60-200	4500X5000X5000
YS-FD-400BG	1300	100-400	5000X5000X6000
YS-FD-600BG	2000	200-600	6000X6000X7000

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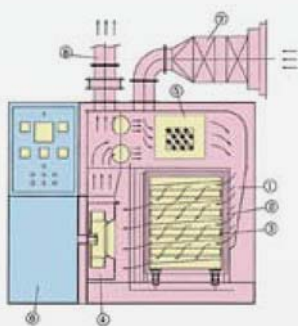
### 流程圖 | FLOWCHART +

1. Drying Chamber
2. Drying Trolley
3. Drying Tray
4. Circulation Fan
5. Heat Exchanger
6. Control Panel
7. Fresh Air Inlet
8. Exhaust Damper

### THEORY +

The drying product is placed on the drying tray which are made of special screen or punched hole plate. The drying tray is placed in the active trolley which has several layers, hot air pierces into the product vertically and removes moisture by special guide flow design to get the drying purpose.

### APPLICATION +

Drying for pharmaceutical, Chinese herbal medicine, Health food, any kind of chemical wet granule and powder

### FUNCTION +

Drying for powder, granule, fine granule, long granule, pellet which contains moisture until get required content of moisture.

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型號 Model	實際操作量 (KG) 依假比重0.4計算	外型尺寸 Size (MM) (W/L/H)
YS-TD-10	60-100	900X2200X1950
YS-TD-15	75-150	1200X3000X2250
YS-TD-20	100-200	1400X3200X2500
YS-TD-30	150-300	2200X3500X2500
YS-TD-40	200-400	2700X3500X2800

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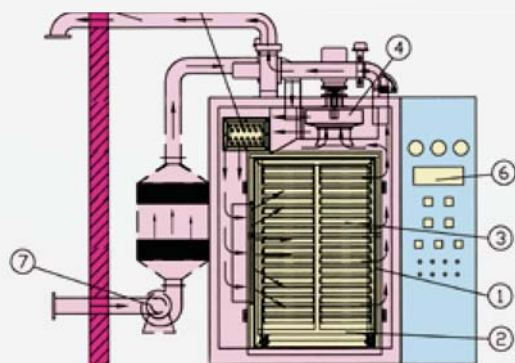
### 流程圖 | FLOWCHART +

- |                    |                    |
|--------------------|--------------------|
| 1. Drying Chamber  | 5. Heat Exchanger  |
| 2. Drying Trolley  | 6. Control Panel   |
| 3. Drying Tray     | 7. Fresh Air Inlet |
| 4. Circulation Fan | 8. Exhaust Damper  |

### THEORY +

The drying product is placed on the drying tray which are made of special screen or punched hole plate. The drying tray is placed in the active trolley which has several layers, hot air pierces into the product vertically and removes moisture by special guide flow design to get the drying purpose.

### APPLICATION +

Drying for pharmaceutical, Chinese herbal medicine, bio-tech product, Health food, any kind of chemical wet granule and powder.

### FUNCTION +

Drying for powder, granule, fine granule, long granule, pellet which contains moisture until get required content of moisture.

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型號 Model	實際操作量 (KG) 依假比重0.4計算	外型尺寸 Size (MM) (W/L/H)
YS-DO-1	60-100	800X2200X1950
YS-DO-2	75-150	800X3000X1950
YS-DO-3	100-200	800X3200X2200
YS-DO-4	150-300	1200X3500X2500
YS-DO-5	200-400	1200X3500X2800

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### THEORY +

Use high speed rotating blades to crush or grind the raw materials to desire size which is smaller than the punch hole, then extrude the granule toward outside of screen, to get the crushing or granulating purpose.

### FUNCTION +

Western medicine, chinese medicine, health food, bio-tech product.

### 規格表 | SPECIFICATION

型號 Model	產量 Capacity (KG)	外型尺寸 Size (MM) (W/L/H)
YS-PM-1	02-05	600X700X1300
YS-PM-3	40-80	800X900X1500
YS-PM-5	50-120	800X900X1600
YS-PM-10	80-200	600X2100X1650
YS-PM-15	150-300	700X2200X1700
YS-PM-20	200-400	800X2400X1800

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## DOUBLE -SHAFT DRY GRANULATOR

+

### THEORY +

Use two closed dry granulate claw to extrude the bigger and irregular granule toward outside of screen, then to form the required even granule.

### FEATURE +

- Multi-function for milling, wet & dry granulation.
- Simple structure is easy to assemble and clean.
- Easy for screen renewing and operating.
- Different adjustable speed for different requirement of product.
- Less dust, low temperature, low noise and easy cleaning.
- Can be integrated with vacuum suction device.
- Special closed shaft design to prevent the black spot and ensure product quality.
- Safety device to protect the operator.
- 360° screen design enables better and faster milling result, output capacity is bigger than other model.
- All the design and manufacture are qualified cGMP standard and FDA requirement.

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### FUNCTION +

- **Crushing**: recover tablet and capsule, or crush the lump into original powder.
- **Milling**: mill the hard or thick raw materials into powder.
- **Dry granulation**: granulate irregular size of granule to even granule after drying.
- **Wet granulation**: granulate the wet powder into granule.

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### THEORY +

- Using high speed rotating blades to crush or grind raw materials to a desired size which is smaller than the perforated plate screen.
- Speed is adjustable to suit different material.
- Perforated plate screen: mesh size can be selected to meet the processing requirements.

### FUNCTION +

- **Crushing**: crush raw material into small granule or powder.
- **Milling**: process larger and irregular granules into smaller and even size.

### FEATURE +

- Special bearing cooling system : for high-speed operation that prevents the high heat damaging the product .
- Suitable for dry and wet granules.
- The safety device will prevent the operator to get hurts.
- Discharge system : the vacuum suction device will be available to eliminate dust contamination. (Optional)

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### THEORY +

The Z-shaped mixer is improved of cone tank, wide in the middle, narrow and unsymmetrical at both ends. Inside have baffles in the middle, when the mixer rotating the products are rolling from top toward bottom and hitting the middle baffles to get the best and efficient mixing result.

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### THEORY +

V-shaped mixer is made of two column tanks, through rotating the granules and separating into two tanks, then two combine to one again, get the best and efficient mixing result.

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### APPLICATION +

Apply for mixing powder, granule, lubricant additive, binder, essence oil of flavor of Western medicine, Chinese medicine and health food.

### FUNCTION +

**Powder mixing:** Mix different kind of powder materials to required uniformity.

**Granule mixing:**

- Mix the granules with lubricant, binder or essence oil of flavor.
- Mix different kind of granules to required uniformity.

### FEATURE +

- Convenient and precise for cleaning.
- Mixing ratio up to 1 / 10,000.
- With safety railing and interlock device.

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### THEORY +

The eight angles shaped tumbler has baffle on each side. The granules are twisted flowing along the angle of baffles and mixed again by the center shaft baffles to get the best and efficient mixing result.

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### THEORY +

The cone shaped mixer is wide in the middle and narrow at both ends. Inside have baffles in the middle, when the mixer rotating the product is rolling from top toward bottom and hitting the middle baffles to get perfect mixing result.

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### THEORY +

Granules or powder are sucked directly from container into mixing machine through conveyor pipe by vacuum. The free powder are collected by sanitary grade dust collection device. It is a device with dust-free, effort-saving, manpower-saving and non-contamination.

### FUNCTION +

Granules or powder are sucked from container of Fluid bed dryer to granulator. The granules after granulation can be conveyed to mixing machine directly.

### APPLICATION +

Pharmaceutical, food, chemical industry, ceramics, fodder, pigments, dyes, agriculture

### FEATURES +

- Reduse dust
- Easy to operation
- Avoid contamination
- saving manpower

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#### THEORY +

- **Lift**: lifting drums by hydraulic cylinder, can be positioned automatically.
- **Flip**: Can be rotated by reducer, and can be 180° positioned automatically.
- **Machine platform rotation**: Can be 90° rotated manually.

#### FUNCTION +

- For Fluid bed dryer/granulator discharge use.
- For drums lifting and feeding.
- For feeding into tablet pressing machine, capsule filling machine, coating machine, etc.

#### APPLICATION +

Pharmaceutical, cosmetic, agriculture, health food or bio-tech products.

#### FEATURES +

- Let powder, granule, and tablet feeding automatically.
- Operation by wire, easy to use and clean.
- Closed operation, reduce dust.
- Can be connected with cone mill, save manpower.

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換桶式  
Change pan type

## THEORY +

Hot airflow pass through a special designed film, heat the uniform rolling tablets inside, spray solution downward on the rolling tablet through pump and spray gun to get coating and drying purpose.

## FUNCTION +

- Film coating: Spray one thin protection layer on the surface of tablet (used in aqueous or organicsolvent ).
- Enteric coating: Spray one protection layer on the surface of tablet to against stomach acid.
- Sugar coating: Multi-layers of sugar coated to protect the tablet.

## APPLICATION +

Pharmaceutical:

- Western medicine: film coating, enteric coating, sugar coating, control released coating.
- Chinese medicine: film coating.

Food:

- Film coating: agriculture products, crystal sugar health food.
- Sugar coating: chewing gun, yeast candy, health food.

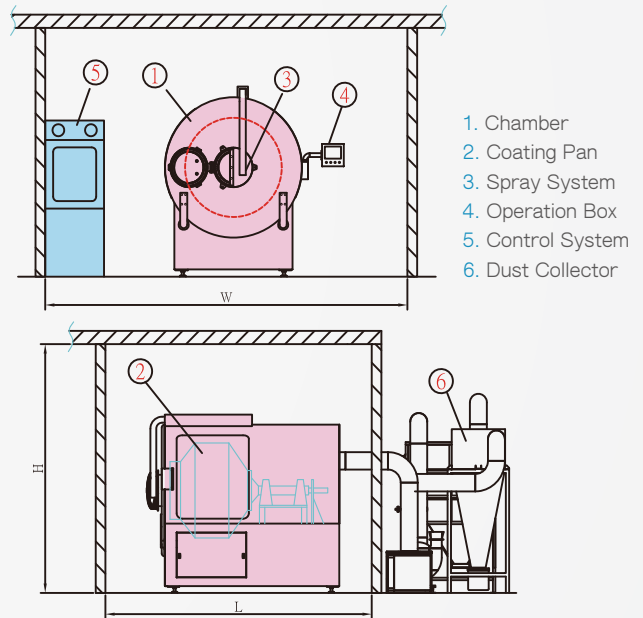
Seed:

- Protection layer coating or add nutrition.

## FEATURE +

- Film and sugar coating are in one machine.
- Punched hole pan design for efficient drying.
- Short operation time:
  - Film coating: 1 ~ 3 hr/batch.
  - Enteric coating: 3 ~ 4 hr/batch.
  - Sugar coating: 6 ~ 16 hr/batch.
- Explosion proof system is available for organic solvent process.
- Auto discharge system for easy discharge.
- Auto control design for sugar coating, even no experienced operator can easy operate also. Well mixed sugar syrup, automatic fixed amount and fixed time, easy operation.
- Uniform coloring, weight and smooth surface for sugar coating.
- Operation time about 6 ~ 16 hr/batch of sugar coating.

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### THEORY +

Place the sugar cores on the turntable. Hot air is blown upward between turntable and granulation area. The air causes the cores to roll, the binder solution is sprayed on the rolling cores through the pump and spray gun, coating and drying simultaneously. If powder coating is required, the fixed volume powder will be sprayed at the same time let solution to coat the powder on the cores, to produce pellet and get coating and drying purpose.

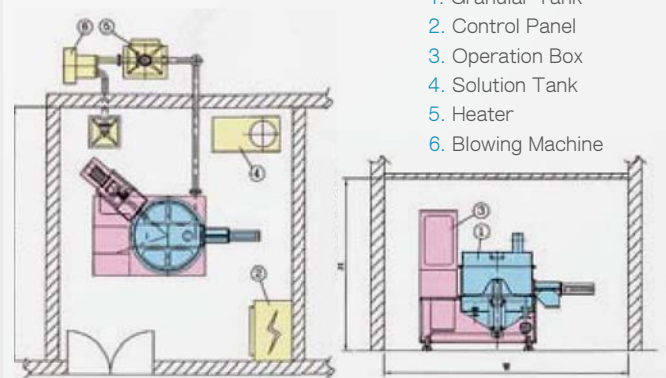
### FUNCTION +

- Powder coating: Binder and powder are sprayed on the tiny granule or sugar to form a bigger pellet.
- Suspension coating: Mixed the active ingredients with binder into solution, spray solution on the smaller pellet to become a bigger pellet.
- Film coating: Spray film / enteric coating on the granulated pellet.

### APPLICATION +

Pharmaceutical, cosmetic, agriculture, health food or bio-tech products.

### 流程圖 | FLOWCHART +

### 規格表 | SPECIFICATION

型號 Model	桶身直徑 Diameter Tank (CM)	實際操作量 Working Batch (KG)	外型尺寸 Size (MM) (W/L/H)
CG-35	35	01-02	2500X2500X2100
CG-75	75	15-30	3000X3000X2300
CG-100	100	30-50	4000X4000X2700
CG-125	125	50-100	5000X5000X3000

INORA

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---

BIO-TECH / CHINESE  
HERBAL MEDICINE  
EQUIPMENT

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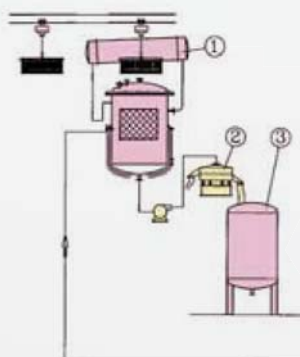
### 流程圖 | FLOWCHART +

1. Extraction Equipment
2. Residue Separator
3. Solution Tank

### THEORY +

The herbs are loaded in the punched hole screen basket and lifted and loaded by crane into the tank, use steam or other heat source to boil, then get into a storage tank after filtering.

The low-temperature concentrator which designed of vacuum theory will evaporate and cool the residual water and alcohol to alcohol water, finally be recovered and used again by distiller.

### APPLICATION +

Extraction of Chinese herbal medicine, Bio-tech product, Health food, which use of water and alcohol.

### FUNCTION +

Use water and alcohol to extract efficient ingredient from natural herbs, then remove and recover residual water and alcohol, preserve the solid materials.

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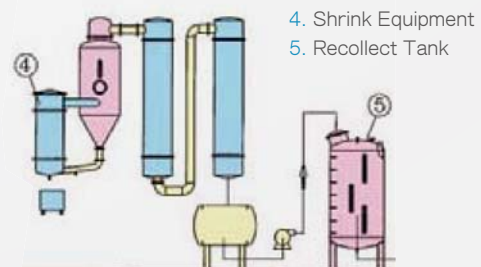
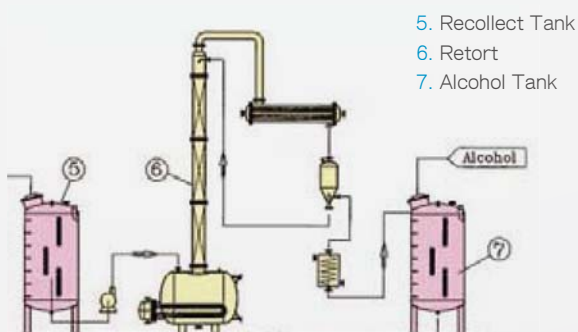
### APPLICATION +

1. This equipment divides into intermission and continuous type. It is consist of distilling kettle, distilling tower body, condenser, buffer tank, flowmeter.
2. Alcohol be recovered and used again by distiller, can reduce cost and protect environment.
3. Continuous type with a pre-heater can be carried out continuous production line with high-volume production needs.
4. It can distill the weak alcohol which purity is lower than 30% purify to 90-95% and the surplus residue alcohol purity 3%.
5. To be equipped with PLC control system and control the parameter of temperature, flux, purity to enhance the production efficiency automatically.

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### APPLICATION +

Chinese medicine, western medicine, bio-tech  
product, health food, food, chemical industry.

### FEATURE +

- Use SUS#304 and 316 stainless steel.
- The size of volume is according to customer requirement.
- The surface contacts directly with product is polished over than 0.1 0.2 Ra.
- All the design and manufacture are qualified cGMP standard and FDA requirement.

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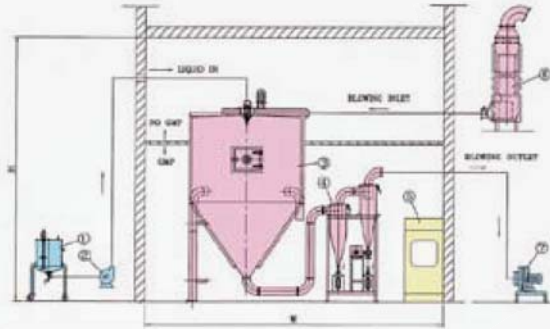
### THEORY +

Heated and purified air flows into the column through exhaust and suction fan, use one set of fast rotating spray nozzle to atomize the solution or paste liquid which sent by pump into very small drop, the drop will combine with heated air and moisture be evaporated to get drying purpose. The original solid ingredient will be collected by collector.

## 流程圖 | FLOWCHART +

- |                   |                    |
|-------------------|--------------------|
| 1. Solution Tank  | 5. Control Panel   |
| 2. Pump           | 6. Heater          |
| 3. Granual Body   | 7. Blowing Machine |
| 4. Dirt Collector |                    |

## APPLICATION +

Chinese herbal medicine, meat seasoning, fruit favored powder, milk powder, chemical raw materials.

- CIP auto cleaning system
- Dehumidifier facility
- All the design and manufacture are qualified cGMP standard and FDA requirement.

## FUNCTION +

Dry into powder from the solution or paste liquid.

## FEATURE +

- Dry into powder from the solution or paste liquid.
- The original essence of material will not be changed after drying.
- Able to dry high viscosity and concentration liquid.
- Instant dry, short drying time.
- Even for hot sensitive materials, also can retain original color, favor and smell.
- Air volume control is proportional by automatic air gate.
- PLC auto control system or touch screen interface for easy operation.
- Explosion proof system is available for organic solvent process.
- Optional equipment for different product feature.
- Temperature recorder

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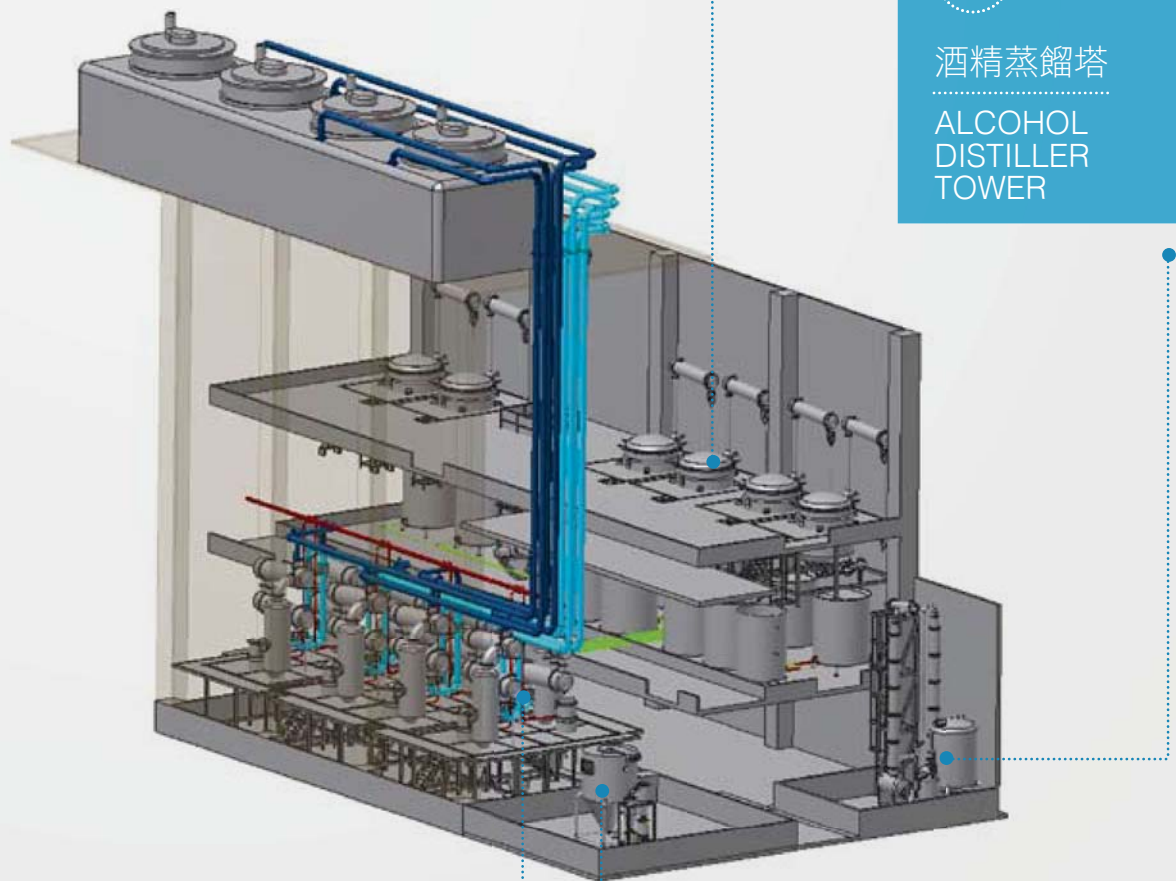
型號 Model	桶身直徑 Diameter of Tank (MM)	產量(L/HR) 依噴霧入水量計算	室內配置尺寸 Size (MM) (W/L/H)
YS-SD-2	800	01-02	2500X3000X2500
YS-SD-5	900	02-05	3000X3500X3000
YS-SD-10	1200	06-10	4000X4000X3500
YS-SD-15	1400	10-15	4500X4500X4000
YS-SD-20	1500	12-20	5000X5000X4500
YS-SD-30	1750	20-30	5000X5500X6000
YS-SD-50	2000	30-50	6000X6000X7000

+

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## PRODUCTION PROCESS FLOW CHART

4

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EXTRACTION  
EQUIPMENT

1

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ALCOHOL  
DISTILLER  
TOWER

3

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VACUUM  
CONCENTRATION  
EQUIPMENT

2

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SPRAY  
DRYER

# INORA

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---

STERILE  
INJECTING  
EQUIPMENT

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## SEMI-AUTOMATIC ROTARY BOTTLE WASHING MACHINE

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### APPLICATION +

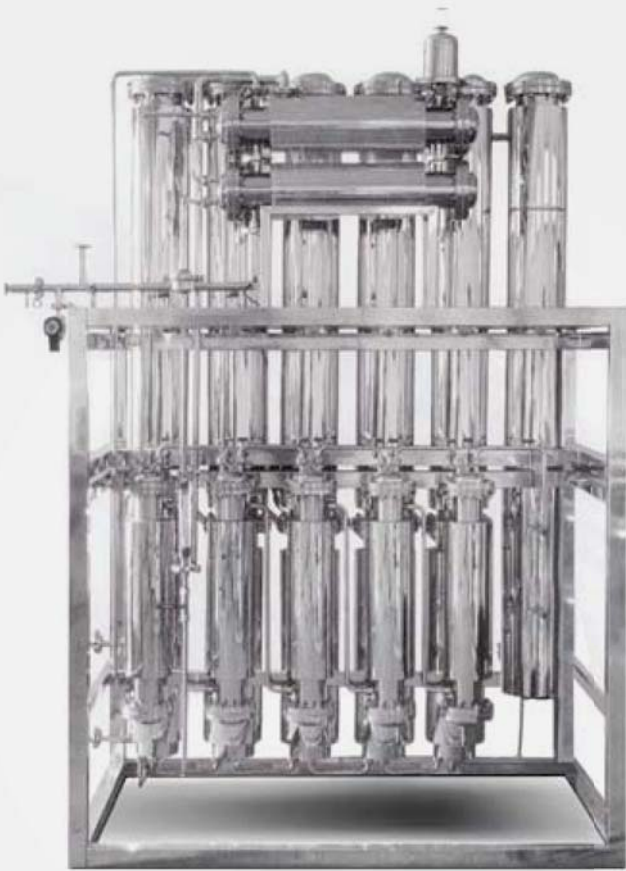
Glass or plastic bottles washing for pharmaceutical、food industry. Washing bottles of various sizes and small capacity

### FUNCTION +

- Equipment is compact, space saving.
- Whole washing process is automatic except manually locate bottles.
- Easy Mould changing to suit various size of bottles.
- Recycling of city water system for energy saving. (Optional)
- Simple operation.
- Adjustable washing speed.

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### FEATURE +

1. Sanitary welding grade: polishing inner and outer for above 0.1 – 0.2 Ra polish.
2. Water/vapor separable design to prevent non-qualified vapor to contaminate the distilled water
3. Distilled water and pure water can be used.
4. Machine contacting feed-water part all using SUS 316L stainless steel ,320-400# grit polish.
5. Double tube sheet design, free from cooling water contamination into distillations by glass fiber.

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6. The distillate temperature is over 80°C to avoid bacteria growing.
7. All designs are according to the world safety pressure vessel standard.
8. This still designed to meet to FDA specification of Water For Injection.

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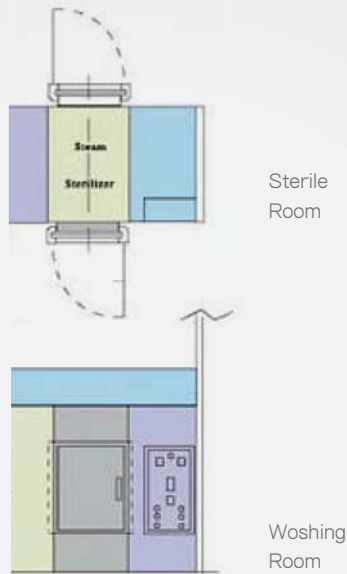
## 流程圖 | FLOWCHART

## APPLICATION +

Pharmaceutical、Food、Hospital、Research Laboratory

## FUNCTION +

The sterilizer is commonly located between the washing room and sterilization room. The extermination of germs and micro-organisms can be achieved when the object for sterilizing is in the contained sterilization room with an essential high temperature and sufficient time; and therefore safer and more reliable products or instruments can be obtained.

## FEATURE +

- Oil hydraulic door: the door clamping device uses an oil hydraulic cylinder to drive clamp for easier and contained operation.
- Steam piping: steam piping fitting is made of sanitary pipe. The steam passes through one micro pre-filter 1  $\mu$  m
- Even steam distribution: by vacuum pump and entrance of steam into sterilization room during sterilization to achieve  $\pm 1$  sterilization temperature.
- Fully automatic control: Adopt PLC control system, when the conditions have been set, the process will be controlled automatically.
- Temperature recorder: records the temperature progress distribution completely and record pressure changes throughout whole sterilization process.
- Inspection hole: validation testing holes are available for inspection.
- Computerized Operation: the computer can calculate the FO figure and shows that there is no more than one germ alive in one million of sterilized objects.
- Three functions in one: integrated sterilization, cooling and leakage test to save time, energy, space and labor.
- Provide complete original record of sensors and calibration of all instruments.

## 規格表 | SPECIFICATION

型號 Model	Volume (L)	Chamber size (WxHxD)	Dimensions (WxLxH)	Steam Consumption (kg/Hr)	Vacuum Pump (HP)	Weight (kg)
YS-SS-041	96	400X400X600	700X1300X900	60	1	1000
YS-SS-061	360	600X600X1000	1000X1300X1300	100	2	1500
YS-SS-081	800	800X1000X1000	1200X1700X1300	130	3	2000
YS-SS-101	1440	1000X1200X1200	1400X1900X1500	200	3	3000
YS-SS-102	2400	1000X1200X2000	1400X1900X2300	280	5	4000
YS-SS-121	1872	1200X1300X1200	1600X2050X1500	240	3	3500
YS-SS-122	3120	1200X1300X2000	1600X2050X2300	350	7 1/2	5000
YS-SS-141	2680	1400X1600X1200	2000X2350X1500	310	5	4500
YS-SS-142	4480	1400X1600X2000	2000X2350X2300	415	10	6000
YS-SS-143	8060	1400X1600X3600	2000X2350X3900	630	15	9000

+

+

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## HOT AIR STERILIZER

### THEORY +

Hot air sterilizer is applicable for sterilizing instruments which can tolerate high temperature and needs to remain in sterilized condition. Under the same temperature, hot air sterilization needs a higher temperature to kill germs and micro-organism than saturated vapor sterilization.

### APPLICATION +

Sterilization of Vial, Ampoule and stainless steel instrument.

### FEATURE +

- Safety sterilized welding: sterilization chamber is welded using stainless steel SUS304 to avoid contamination from the insulation material after long use.
- Complete contained: during the process of sterilization and when the sterilizer is off, the dampers are always tightly closed to ensure the sterilizing chamber is in a contained condition.
- High efficiency filter : the absolute Heap Filter 99.97%, this filter can tolerate temperature above 350°C and filter particles below 0.3 Mirco.

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Washing Room

### 規格表 | SPECIFICATION

型號 Model	Chamber Size (MM) (WXLXH)	Overall Dimensions (KG) (WXLXH)	Maximum Temperature (C)	Ampoule 10CC (P.C.S BATCH)	Vial 10CC (P.C.S BATCH)	Vial 20CC (P.C.S BATCH)	Vial 100CC (P.C.S BATCH)
YS-DHS-1	810X1100X1000	1600X2500X1300	280	10000	11000	5300	1500
YS-DHS-2	1600X1500X800	2600X2880X1100	280	18000	19000	9900	300
YS-DHS-3	1600X1500X1330	2600X2880X1630	280	27000	28000	13000	5000

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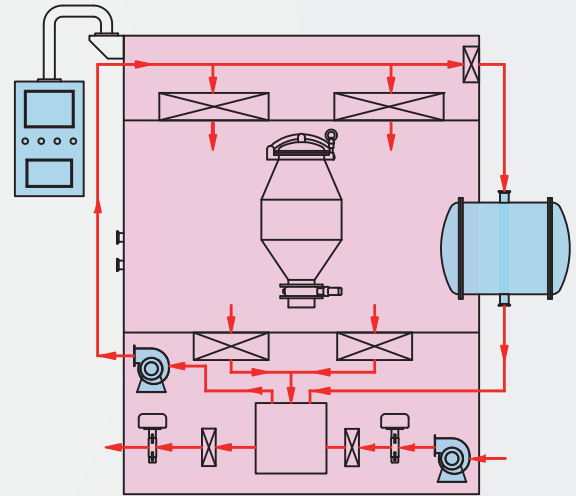
## FEATURE +

- Glove recommended for polluting handling with high mechanical and chemical risks.
- Recommended for pharmaceutical and electronics industries, chemical and medical research.
- High flexibility, dexterity and tactility.
- Excellent chemical resistance to mineral and organic acids, alcohols, oils, lubricants and hydrocarbons.
- Operating temperature range : -30°C to +50°C
- Dipped in 100% neoprene.
- PLC automated control system
- Interlocks on inlet/outlet doors.
- Inlet/ outlet HEPA filters.
- All the design and manufacture are qualified PIC/S cGMP standard and FDA requirement.

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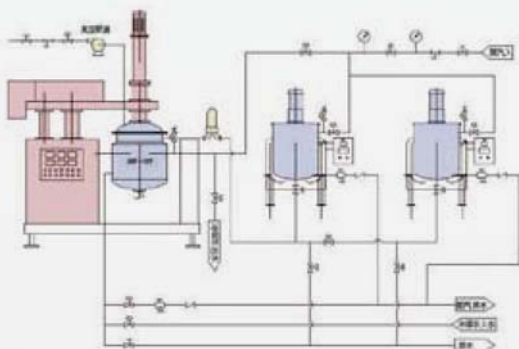
### 流程圖 | FLOWCHART +

### APPLICATION +

Pharmaceutical: ointment.  
Cosmetic: cream, toothpaste, etc.

### FEATURE +

- Water, oil solution mixing tank with heating system for convenient of raw material heating and dissolving.
- Vacuum system: the oil and water solution can be conveyed to emulsifying tank by vacuum transfer in a closed operation.
- Set variables during operation: heating temperature, homogeneous speed, scraper speed.
- Automatic process: heating / emulsify / vacuum / cooling.
- Oil hydraulic system: lift up the scraper with cover for easy cleaning.
- To use operator interface ( process visualization ) control system for the whole process automation.

### 規格表 | SPECIFICATION

型號 Model	全容積 Full Volume (L)	最大高度 Max. Height (MM)	外型尺寸 Size (MM) (WXLXH)
YS-HM-100	100	3450	750X2000X2450
YS-HM-150	150	3500	800X2100X2500
YS-HM-200	200	3600	870X2350X2600
YS-HM-250	250	3700	900X2400X2650
YS-HM-350	350	3800	930X2440X2700
YS-HM-500	500	4350	1350X2800X3000